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INSTRUCTION MANUAL

ZOOM STEREO MICROSCOPE SMZ-0790

Thank you for purchasing the Zoom Stereo Microscope SMZ-0790. Please read this instruction manual thoroughly before use to ensure proper operation and long service.


NOTES

- This is a precision instrument – Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Avoid direct sunlight or areas with high temperatures or high levels of dust.
Keep in location where temperature and humidity are stable.
- Keep lens surface clean from dirt and fingerprints.
- Clean the lens gently with lens cleaning cloth or soft cotton cloth.
- Do not turn the left and right Zoom Controls in opposite directions; this will damage the instrument.
- When not in use store with protective cover, in location that has low moisture in order to prevent internal mold growth.

SPECIFICATIONS

Magnification	7~45× • 14~90×
Objective Lens	0.7~4.5× (Zoom Ratio 1:6.4)
Eyepiece Lens	WF10× (20mm Field of View) WF20× (10mm Field of View)
Working Distance	100mm
Eyepiece Angle	45°
Focus Adjustment	Movable Body
Stand	Pole Type (Φ32mm)
Stage Plate	Φ95mm Black/White Plastic Plate
Eye Width Adj. Range	55~75mm
Diopter Adj. Range	±5 Diopter
Illumination	Optional Ring Light (sold separately)
Base Dimensions	L275 × W205 × H22mm
Weight	approx. 4.0kg

PART IDENTIFICATION and ASSEMBLY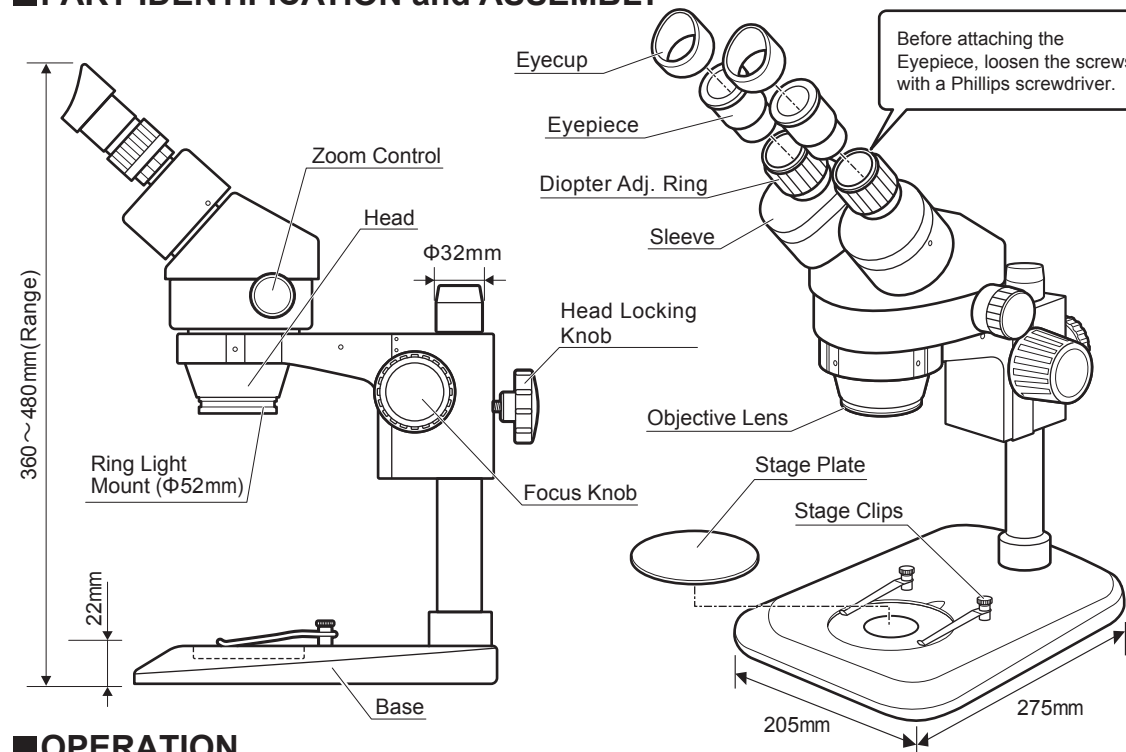

OPERATION

1. Sample Placement

- ① Place the sample in the center of the Stage Plate.
- ② If required, use Stage Clips to hold sample in place.

2. Eyecup Use

- **For use without eyeglasses:**
Hold the Diopter Adj. Ring to prevent it from rotating, turn the eyepiece, align the eye cup with the observer's face, and bring the eye close to observe.

- **For use with eyeglasses:**
Remove Eyecups to use.

3. Eye Width Adjustment

Hold the left and right sleeve in each hand and look through the Eyepieces. Adjust the Sleeves to align the left and right field of view seen through the Eyepieces.

4. Focus Adjustment•Diopter Adjustment

- ① Rotate the Zoom Control to set for highest magnification.
- ② With right eye, look through the right eyepiece and adjust the focus using the Focus Knob.

- ③ Rotate the Zoom Control for lowest magnification.
- ④ Using the Right Diopter Adj. Ring, adjust the focus if required.
- ⑤ Now rotate the Zoom Control back to highest magnification and check the focus.
Repeat steps ② to ④ to achieve the best focus accuracy.
- ⑥ Look through the left Eyepiece with the left eye and adjust focus using the Left Diopter Adj. Ring similar to steps ① to ⑤.

5. Focus Knob Tension Adjustment

If the Focus Knob movement is too stiff or too loose, adjust the knob tension by holding either side of the Focus Knobs in place and turning the other side.
Turn it clockwise to add the knob tension, then eventually the head position is fixed. Turn it counterclockwise to reduce the knob tension.

Niigata seiki Co., Ltd.

5-3-14, Tsukanome, Sanjo, Niigata, Japan, 955-0055
Tel. : +81-256-33-5522 Fax. : +81-256-33-5518
MAIL intl.sales@niigataseiki.co.jp
URL http://www.niigataseiki.co.jp